

## Initial set-up for Scanning Tunnelling Microscope



In order to avoid driver conflicts, it is important to follow the following steps carefully.

Connect the USB cable to the control unit but do not yet plug it into the computer.

Connect the control unit to its power supply but do not turn it on yet.

Set up the scan head and connect it to the control unit.

Switch on the computer and start Windows.

Insert the install CD and run the software installation.

Click the button that says "Install Easy Scan 2 Software" and go through the installation procedure to completion.

Switch on the control unit and connect to the computer via the USB cable.

Wait for the drivers to install.