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U14600 Optics on magnetic boards, basic set

Instruction sheet

3/11 ALF

The demonstration equipment set (basic set) is for the conducting of numerous demonstration experiments in geometric optics, e.g. casting of shadows, laws of reflection, laws of refraction, total reflection, angle of least deflection in prisms, focal lengths of mirrors and lenses, laws of lenses and imaging errors.

1. Safety instructions

- Beware. Lights become hot when left on for a long period.
- Do not clean optical elements with highly active fluids or solvents. This may result in damage.

2. Description, technical data

All optical elements in the basic set are coated with magnetic foil and are available separately.

Height: 35 mm each

Plane mirror

Metal strip on plastic block
Dimensions: 200 mm x 35 mm

Convex and concave mirrors

2 metal strips on plastic block
Dimensions: 200 mm x 35 mm
Radius of curvature: 200 mm
Focal length: ± 100 mm

Plano-concave lens

Transparent acrylic body with matt surface
Dimensions: 200 mm x 40 mm x 35 mm
Radius of curvature: 200 mm
Focal length: -400 mm

Plano-convex lens

Transparent acrylic body with matt surface
Dimensions: 200 mm x 40 mm x 35 mm
Radius of curvature: 200 mm
Focal length: 400 mm

Semi-circular lens

Transparent acrylic body with matt surface
Radius of curvature: 100 mm
Focal length: 200 mm

Plane parallel block

Transparent acrylic body with matt surface
Dimensions: 200 mm x 100 mm x 35 mm

Right-angled prism

Transparent acrylic body with matt surface
Dimensions: 200 mm x 200 mm x 283 mm

Set of three opaque bodies for casting shadows

Plastic forms
Cuboid: 100mm x 20 mm x 35 mm
Cylinder: 60 mm \varnothing x 35 mm
Cylinder: 15 mm \varnothing x 35 mm

2.1 Scope of delivery:

No.	Cat. no.	Description
1	U15510	Plane mirror, magnetic
1	U15511	Convex and concave mirrors, magnetic
1	U15515	Plano-concave lens, magnetic
1	U15516	Plano-convex lens, magnetic
1	U15517	Plane-parallel block, magnetic
1	U15518	Semi-circular lens, magnetic
1	U15520	Right-angled prism, magnetic
1	U15525	Set of three opaque bodies, magnetic

2.2 Accessories:

- 1 Magnetic board or whiteboard (U10030 or U10031)
 - 2 Single-ray projectors U40120
 - 2 Magnetic holders for single-ray projector U40121 or
 - 1 Multiple ray projector U40110
- Power supply, 12 V, 10 A, e.g. U11735

3. Cleaning

- Clean optical elements using a damp cloth with washing-up liquid or plexiglass cleaner.

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